



Biogal's **ImmunoComb[®]**

Mouse and Murine Antibody Test Kits

Convenient Solid Phase ELISA Antibody Test Kits.

The kits are completely self-contained.

They require no specialized space, equipment or personnel.

The ImmunoComb[®] Mouse and Murine kits were developed through collaborative research with Charles River Laboratories (CRL). CRL thought that the ImmunoComb[®] technology could be applied for in-house health monitoring of laboratory animal colonies. They wanted to develop a kit for facilities that did not have fully equipped diagnostic laboratories. The Murine and Mouse ImmunoComb[®] Antibody Test Kits was the outcome of this project.





The ImmunoComb® Mouse and Murine Kits offered by Biogal Galed Laboratories are convenient solid phase ELISA Antibody Test Kits.

- ✓ The kits are completely self-contained.
- ✓ They require no specialized space, equipment or personnel.

The kits detect antibodies to:

- 1. Sendai Virus, Rodent Coronaviruses and Mycoplasma pulmonis (Murine Kit)**
- 2. Mouse Hepatitis Virus, Mouse Parvovirus (strains 1 and 2), Minute Virus of Mouse and Murine Norovirus (Mouse kit).**

Advantages:

- ✓ Reagents need no preparation
- ✓ Plain tap water is used for wash steps
- ✓ No incubator is needed
- ✓ Results are ready in 90 minutes (hands-on time is much less)
- ✓ Economical
- ✓ Low cost per assay
- ✓ Each kit supplies 3 combs and 3 plates for 90 determinations on 30 samples
- ✓ Comb breaks apart to test as few as one sample

Life Technologies (India) Pvt. Ltd.

Tel # 91-11-4220-8000; 4220-8111; 4220-8222, Fax# 91-11-4220-8444, Mobile# +91-98105-21400,
Email - customerservice@lifetechindia.com Web - www.lifetechindia.com

ImmunoComb® Mouse and Murine Kit Contents:

Both of the ImmunoComb® Kits are completely inclusive and they do not require specialized space, equipment or training. The reagents included need no preparation and only water is used for the wash steps.



Kit contents (pictured above) include:

1. ImmunoComb® Cards, 2. Developing Plates, 3. CombScale, 4. Control Sera, 5. Heparinized Capillary Tubes (5a) and Wire Piston (5b) – additional cost, 6. Tweezers, 7. Instructions for Use and Results Sheet (not pictured)

The "CombScan" Software:

The "CombScan" Software is a computer program that can read the ImmunoComb results faster and in a more accurate manner. This is a free of charge option for customers that need it. Biogal provide the disc and calibration comb and customers need a computer and appropriate scanner.


To download the complete ImmunoComb Mouse and Murine Product Information and Instruction Manual, Please visit the following link:

http://www.lifetechindia.com/Life_Technologies_India_ImmunoComb_Mouse_Murine_Kits.html


Life Technologies (India) Pvt. Ltd.

Tel # 91-11-4220-8000; 4220-8111; 4220-8222, Fax# 91-11-4220-8444, Mobile# +91-98105-21400,
Email - customerservice@lifetechindia.com Web - www.lifetechindia.com

Directions For Use:

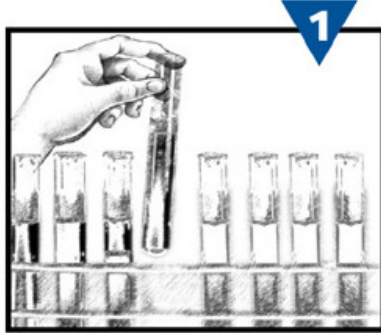


1) Obtain blood/serum from animal/s

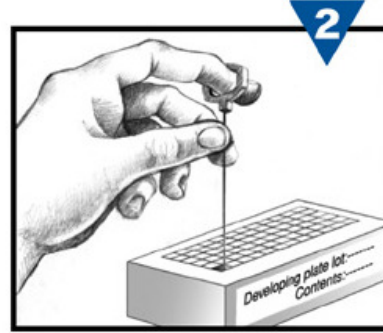


2) Deposit sample/s into well/s A in developing plate



3) Insert Comb into well/s A. Incubate for a specified time

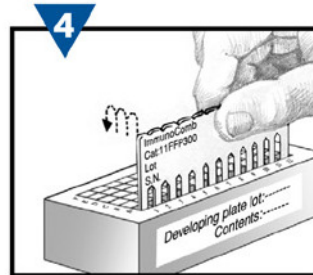

4) Transfer the Comb from well/s A to well/s F in a timed sequence, according to specific kit.

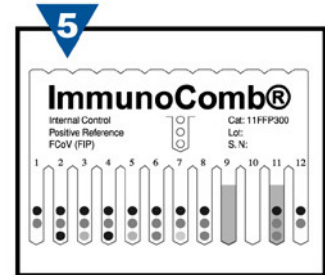

5) Let comb dry and Read results. The color intensity of spots on the comb corresponds to antibody level.

Life Technologies (India) Pvt. Ltd.

Tel # 91-11-4220-8000; 4220-8111; 4220-8222, Fax# 91-11-4220-8444, Mobile# +91-98105-21400,
Email - customerservice@lifetechindia.com Web - www.lifetechindia.com