

## MG1655 Chemically Competent Cells

Catalog #	Package Size
1076-12	12x50 µl

### Description

Intact Genomics (ig®) chemically competent MG1655 chemically competent cells are suitable for routine DNA transformation and other research purposes. MG1655 is the "wild type" K-12 strain with minimal genetic manipulation having only been cured of the temperate bacteriophage lambda and F plasmid by ultraviolet light and acridine orange (1).

### Specifications

Competent cell type: Chemically competent  
Derivative of: MG1655  
Species: *E. coli*  
Format: Tubes  
Transformation efficiency:  $\geq 5.0 \times 10^6$  cfu/µg pUC19 DNA  
Blue/white screening: No  
Shipping condition: Dry ice

### Reagents Needed for One Reaction

ig® MG1655 Chemically Competent Cells: 50 µl  
DNA (or pUC19 Control, 10 pg/µl): 1 µl  
Recovery medium: 1 ml

### Product Includes & Storage

- ig® MG1655 Chemically Competent Cells: -80 °C
- pUC19 control DNA: -20 °C
- Recovery medium: 4 °C

### Genomic Features

ig® MG1655 chemically competent cells have the following genomic features:

- Cured of the temperate bacteriophage lambda and F plasmid
- A frameshift at the end of rph
- ilvG- mutation is a frameshift that knocks out acetohydroxy acid synthase II
- rfb-50 mutation is an IS5 insertion that results in the absence of O-antigen synthesis

### Genotype

F- lambda- *ilvG- rfb-50 rph-1*

### General Guidelines

Follow these guidelines when using MG1655 chemically competent *E. coli*.

- Handle competent cells gently as they are highly sensitive to changes in temperature or mechanical lysis caused by pipetting.
- Thaw competent cells on ice, and transform cells immediately following thawing. After adding DNA, mix by tapping the tube gently. Do not mix cells by pipetting or vortexing.

### Example Calculation of TE

Transformation Efficiency (TE) is defined as the number of colony forming units (cfu) produced by transforming 1 µg of plasmid into a given volume of competent cells.

$TE = \text{Colonies}/\mu\text{g}/\text{Dilution}$

Transform 1 µl of (10 pg/µl) pUC19 control plasmid into 50 µl of cells, add 950 µl of Recovery Medium. Dilute 10 µl of this in 990 µl of Recovery Medium and plate 50 µl.

Count the colonies on the plate the next day. If you count 100 colonies, the TE is calculated as follows:

$$\begin{aligned} \text{Colonies} &= 100 \\ \mu\text{g of DNA} &= 0.00001 \\ \text{Dilution} &= 50/1000 \times 10/1000 = 0.0005 \\ \text{TE} &= 100/.00001/.0005 = 2.0 \times 10^{10} \end{aligned}$$

### High Efficiency Transformation Protocol

Use this procedure to transform ig® MG1655 chemically competent cells. We recommend verifying the transformation efficiency of the cells using the pUC19 control DNA supplied with the kit. Do not use these cells for electroporation.

- Remove competent cells from the -80 °C freezer and thaw completely on wet ice (10-15 minutes).
- Aliquot 1-5 µl (1 pg-100 ng) of DNA to the chilled microcentrifuge tubes on ice.
- When the cells are thawed, add 50 µl of cells to each DNA tube on ice and mix gently by tapping 4-5 times. For the pUC19 control, add 1 µl of (10 pg/µl) DNA to a chilled microcentrifuge tube, prior to adding 50 µl of cells. Mix well by tapping. Do not pipette up and down or vortex to mix, this can harm cells and decrease transformation efficiency.
- Incubate the cells with DNA on ice for 30 minutes.
- After 30 minute ice incubation, heat shock the cells at 42 °C for 45 seconds.
- Transfer the tubes to ice for 2 minutes.
- Add 950 µl of Recovery Medium or any other medium of choice to each tube.
- Incubate tubes at 37 °C for 1 hour at 210 rpm.
- Spread 50 µl to 200 µl from each transformation on Pre-warmed selection plates. We recommend plating two different volumes to ensure that at least one plate will have well-spaced colonies. For the pUC19 control, plate 50 µl on an LB plate containing 100 µg/ml ampicillin. Use sterilized spreader or autoclaved ColiRoller™ plating beads to spread evenly.
- Incubate the plates overnight at 37 °C.

## MG1655 Chemically Competent Cells



### 5 Minute Transformation Protocol

The following procedure results in only ~10% of the transformation efficiency as the protocol listed above.

- 1) Remove competent cells from the -80 °C freezer and thaw in your hand.
- 2) Aliquot 1-5 µl (1 pg-100 ng) of DNA to the microcentrifuge tubes. **Do not** pipette up and down or vortex to mix, this can harm cells and decrease transformation efficiency.
- 3) Incubate the cells with DNA on ice for 2 minutes.
- 4) After 2 minute ice incubation, heat shock the cells at 42 °C for 45 seconds.
- 5) Transfer the tubes to ice for 2 minutes.
- 6) Add 950 µl of Recovery Medium at room temperature or any other medium of choice to each tube. Immediately spread 50 µl to 200 µl from each transformation on pre-warmed selection plates. We recommend plating two different volumes to ensure that at least one plate will have well-spaced colonies. For the pUC19 control, plate 50 µl on an LB plate containing 100 µg/ml ampicillin. Use sterilized spreader or autoclaved ColiRoller™ plating beads to spread evenly.
- 7) Incubate the plates overnight at 37 °C.

### Related Products

- ig® 5-Alpha Chemically Comp. Cells (Cat.# 1031-12)
- DirectPlate™ DH5-Alpha Chem. Comp. Cells (Cat# 1013-12)
- Quick10™ Cloning Kit (Cat#4122)
- T4 DNA Ligase (Cat.# 3212)
- i7® High Fidelity DNA Polymerase (Cat.# 3254)

### Technical Support

Intact Genomics is committed to supporting the worldwide scientific research community by supplying the highest quality reagents. Each new lot of our products is tested to ensure they meet the quality standards and specifications designated for the product. Please follow the instructions carefully and contact us if additional assistance is needed. We appreciate your business and your feedback regarding the performance of our products in your applications.

### Quality Control

Transformation efficiency is tested by using the pUC19 control DNA supplied with the kit and the high efficiency transformation protocol listed below. Transformation efficiency should be  $\geq 5 \times 10^6$  CFU/µg pUC19 DNA. Untransformed cells are tested for appropriate antibiotic sensitivity.

### References:

1. Blattner, et al (1997) The complete genome sequence of *Escherichia coli* K-12. *Science* 277(5331), 1453-1462.
2. Jensen (1993) The *Escherichia coli* K-12 "wild types" W3110 and MG1655 have an *rph* frameshift mutation that leads to pyrimidine starvation due to low *pyrE* expression levels. *J Bacteriol* 175(11), 3401-3407.
3. Lawther, et al. (1982) DNA sequence fine-structure analysis of *ilvG* (*IlvG*+) mutations of *Escherichia coli* K-12. *J Bacteriol* 149(1), 294-298.
4. Liu and Reeves (1994) *Escherichia coli* K12 regains its O antigen. *Microbiology* 140(Pt 1), 49-57.
5. Bachmann (1996) Derivations and Genotypes of Some Mutant Derivatives of *Escherichia coli* K-12. p2460-2488 in Neidhardt (ed), *Escherichia coli* and *Salmonella*, 2nd Edition ASM Press, Washington DC

India Contact:  
Life Technologies (India) Pvt. Ltd.  
Ph: +91-11-42208000, 42208111, 42208222  
Mobile: +91-9810521400  
Email: [customerservice@lifetechindia.com](mailto:customerservice@lifetechindia.com)  
[www.lifetechindia.com](http://www.lifetechindia.com)