


Importance of STEC Determination

Shiga toxin-producing *E. coli* (STEC) are food-borne pathogens that cause gastrointestinal illness, hemorrhagic colitis, and hemolytic uremic syndrome. As few as 100 cells can cause disease leading to serious illness or death. *Escherichia coli* O157:H7 is the STEC serotype most often implicated in outbreaks; however, there are numerous other STEC serotypes that have caused serious human illness and outbreaks. STEC that cause human illness were added as notifiable pathogens to the Nationally Notifiable Diseases Surveillance System in 2000. From 2000-2010, 7695 cases were reported; 5688 were associated with the serogroup O157, and the 83% of the other STEC were serogroups O26, O45, O103, O111, O121 and O145. However, other serogroups, including O91, O113, O104, have also caused serious human illness. *Escherichia coli* from each of the serogroups listed above have been found in ground beef, cattle hides, and in feces. Bovine feces can be a source of environmental contamination (soil, water) which can lead to secondary contamination of produce growing in fields.

It is difficult to distinguish pathogenic *E. coli* strains from non-pathogenic *E. coli* strains because the former rarely possess any distinguishing phenotypic or biochemical characteristics which differ from the latter. Methods such as the Shiga toxin ELISA described in this User's Guide have been developed to determine the presence of virulence markers, and therefore the presence of pathogenic *E. coli*. Rapid, sensitive detection that can distinguish between Stx 2 and Stx 1 toxin types are vital for proper diagnosis.

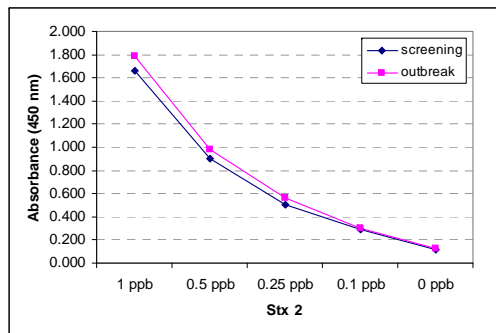

Example of absorbances generated by different concentrations of Stx 2 in media incubated according to section D. **Not intended for use in sample evaluation.**

References

USDA-FSIS Microbiological Laboratory Guidebook Chapter 5.08 "Detection, Isolation and Identification of *Escherichia coli* O157:H7 from Meat Products and Carcass and Environmental Sponges"

General Limited Warranty: Abraxis LLC warrants the products manufactured by the Company, against defects and workmanship when used in accordance with the applicable instructions for a period not to extend beyond the product's printed expiration date. **Abraxis makes no other warranty, expressed or implied. There is no warranty of merchantability or fitness for a particular purpose.**

For ordering or technical assistance contact: Abraxis LLC
54 Steamwhistle Drive
Warminster, PA 18974
Tel.: (215) 357-3911
Fax: (215) 357-5232
Email: info@abraxiskits.com
WEB: www.abraxiskits.com

R0102615

Shiga toxin 2 ELISA (Microtiter Plate)

Enzyme-Linked Immunosorbent Assay for the Determination of Shiga toxin 2 (Stx 2) in Ground Beef Samples



Product No. 542010

1. General Description

The Abraxis Shiga toxin 2 ELISA is an immunoassay for the qualitative screening of Shiga toxin 2 (Stx 2). This test is suitable for the qualitative screening of Stx 2 in samples (please refer to the appropriate technical bulletins for additional matrices procedures). Shiga toxin is a virulence marker for O157 and non-O157 Shiga toxin producing *Escherichia coli* (STEC). This ELISA allows for the rapid and sensitive detection of Stx 2 and when used in combination with Stx 1 ELISA allows the differentiation of the two toxin types.

2. Safety Instructions

The positive control in the test kit contains small amounts of Shiga toxin 2. In addition, the substrate solution contains tetramethylbenzidine and the stop solution contains diluted sulfuric acid. Avoid contact of stopping solution with skin and mucous membranes. If these reagents come in contact with skin, wash with water. Excess beef and incubated media should be considered biological waste, and decontaminated, along with any contaminated equipment, by autoclaving or another effective method.

3. Storage and Stability

The Stx 2 ELISA should be stored in the refrigerator (4–8°C). The solutions must be allowed to reach room temperature (20–25°C) before use. Reagents may be used until the expiration date on the box.

4. Test Principle

The test is an indirect ELISA based on the recognition of Shiga toxin 2 by specific antibodies. Ground beef samples are incubated with a medium that encourages *E. coli* growth, if present, and shiga toxin production. If Stx 2 is present in a sample, is bound by an antibody immobilized on the wells of the microtiter plate. After a washing step, a second antibody is added which binds to the Stx 2 bound to the wells. A second washing step is followed by the addition of an HRP labeled antibody which binds to the existing antigen/antibody complex in the wells. After a final wash and addition of the substrate solution, a color signal is generated. The intensity of the blue color is proportional to the concentration of Stx 2 present in the sample. The color reaction is stopped after a specified time and the color is evaluated using an ELISA reader. The result of the evaluation is compared to a known value to determine whether the sample is positive or below the limit of Stx 2 detection.

5. Limitations of the Stx 2 ELISA, Possible Test Interference

Although many organic and inorganic compounds commonly found in samples have been tested and found not to interfere with this test, due to the high variability of compounds that might be found in samples, test interferences caused by matrix effects cannot be completely excluded.

Samples must be diluted and incubated as instructed in the sample preparation section (Section D) or appropriate technical bulletin before testing in the ELISA.

Mistakes in handling the test also can cause errors. Possible sources for such errors include: improper storage of the test kit, incorrect pipetting sequence or inaccurate volumes of the reagents, too long or too short incubation times during the immune and/or substrate reactions, extreme temperatures (lower than 10°C or higher than 30°C) during the test performance.

The Abraxis Stx 2 ELISA kit provides screening results only and does not differentiate between shiga toxin producing bacteria. Any positive results should be confirmed by PCR or another approved method.

A. Reagents and Materials Provided

1. Microtiter plate (8 wells X 12 strips) coated with a capture antibody, in a resealable aluminum pouch
2. Stx 2 positive control as Stx 2 (1): 0.5 ng/mL (ppb), X mL
3. Stx 2 negative control, X mL
4. Anti-Stx 2 Antibody Solution, X mL
5. Secondary antibody HRP Conjugate, X mL
6. Wash Solution (5X) Concentrate, 100 mL, must be diluted before use, see Test Preparation (Section C)
7. Substrate (Color) Solution (TMB), X mL
8. Stop Solution, X mL (handle with care)

B. Additional Materials (not delivered with the test kit)

1. Micro-pipettes with disposable sterile plastic tips (20-1000 µL)
2. Multi-channel pipette (50-250 µL) or stepper pipette with disposable sterile plastic tips (50-250 µL)
3. Microtiter plate reader (wave length 450 nm)
4. Sterile clear polypropylene bags with mesh (ca. 24" x 30-36")
5. Stomacher (optional)
6. Modified tryptone soya broth with Novobiocin (mTSB+n) or other approved sample incubation medium
7. *E. coli* positive control (for inoculating media)
8. Incubator, static 42 ± 1°C
9. Automatic pipettor with 25 mL sterile disposable serological pipets
10. Paper towels or equivalent absorbent material
11. Timer
12. Autoclave
13. 50 mL sterile disposable conical vials

C. Test Preparation

Micro-pipetting equipment and pipette tips for pipetting the standards and the samples are necessary. In order to equalize the incubation periods on the entire microtiter plate, a multi-channel pipette or a stepping pipette is recommended for adding the enzyme conjugate, antibody, substrate, and stop solutions. Please only use the reagents and standards from one package lot in one test, as they have been adjusted in combination.

1. Adjust the microtiter plate and the reagents to room temperature before use.
2. Remove the number of microtiter plate strips required from the aluminum pouch. The remaining strips are stored in the aluminum pouch and zip-locked closed. Store the remaining kit in the refrigerator (4-8°C).
3. The standard solutions, antibody, conjugate, substrate and stop solutions are ready to use and do not require any further dilutions.
4. Dilute the Wash Solution (5X) Concentrate at a ratio of 1:5. If using the entire bottle (100 mL), add to 400 mL of deionized or distilled water and mix thoroughly.
5. The stop solution must be handled with care as it contains diluted H₂SO₄.

D. Sample Preparation

Beef Samples

(see also USDA FSIS MLG Chapter 5.08: Detection, Isolation and Identification of *Escherichia coli* O157:H7 from Meat Products and Carcass and Environmental Sponges)

NOTE: The Abraxis Stx 2 ELISA may be used with either beef screening samples (sample diluted 1:4 in media) OR outbreak related samples (sample diluted 1:10 in media)

1. Disinfect surface of sample packaging prior to opening
2. i. **Screening samples:** Prepare in a sterile strainer bag a 1:4 dilution (one part beef in three parts media, e.g. 25 g beef with 75 mL enrichment medium). Increase or decrease total volume incubated according to amount needed to prepare representative samples and perform all analyses.
ii. **Outbreak related samples:** Prepare in a sterile strainer bag a 1:10 dilution (one part beef in 9 parts media, e.g. 10 g beef with 90 mL enrichment medium). Increase or decrease total volume incubated according to amount needed to prepare representative samples and perform all analyses.
3. Pummel or hand massage until clumps are dispersed.
4. The USDA requires that both a known positive sample and an uninoculated medium control are incubated alongside each group of samples.
5. Incubate all bags for 15-22 hr at 42 ± 1°C.
6. After removing samples from incubator, mix gently by hand massaging to ensure uniformity and draw off 10-20 mL media from opposite side of mesh from beef sample.
7. Allow to settle 10-15 minutes in 50 mL conical vial. Gently decant top layer to a clean vial if desired, use as sample for ELISA. No further dilution is necessary.

E. Working Scheme

The microtiter plate consists of 12 strips of 8 wells, which can be used individually for the test. **The FDA requires the incubation of O157 inoculated and uninoculated media alongside beef samples; positive and negative controls must be run with each test.** Never use the values of controls that have been determined in a test performed previously.

(+) Ctrl, (-) Ctrl: kit controls

EC (+), EC (-): incubated positive

and negative controls

Samp 1, Samp 2: incubated samples

	1	2	3	4	5	6	7	8	9	10	11	12
A	(+) Ctrl	Samp 1										
B	(+) Ctrl	Samp 1										
C	(-) Ctrl	Samp 2										
D	(-) Ctrl	Samp 2										
E	EC (+)	etc.										
F	EC (+)	etc.										
G	EC (-)											
H	EC (-)											

F. Assay Procedure

1. Add 100 µL of the **control solutions or samples** (Section D) into the wells of the test strips according to the working scheme given. Analysis in duplicate or triplicate is recommended. Cover the wells with parafilm or tape and mix the contents by moving the strip holder in a circular motion on the benchtop for 60 seconds. Be careful not to spill the contents.
2. Incubate for 60 minutes at room temperature.
3. Remove the covering and decant the contents of the wells into a sink. Wash the strips **three times** using the **diluted washing buffer solution**. Please use at least a volume of 250 µL of washing buffer for each well in each washing step. Remaining buffer in the wells should be removed by patting the inverted plate dry on a stack of paper towels
4. Add 100 µL of **Anti-Stx 2 antibody solution** to the individual wells successively using a multi-channel pipette or a stepping pipette. Cover the wells with parafilm or tape and mix the contents by moving the strip holder in a circular motion on the benchtop for 60 seconds. Be careful not to spill the contents.
5. Incubate 30 minutes at room temperature.
6. Remove the covering and decant the contents of the wells into a sink. Wash the strips **three times** using the **diluted washing buffer solution**. Please use at least a volume of 250 µL of washing buffer for each well in each washing step. Remaining buffer in the wells should be removed by patting the inverted plate dry on a stack of paper towels
7. Add 100 µL of **enzyme conjugate solution** to the individual wells successively using a multi-channel pipette or a stepping pipette. Cover the wells with parafilm or tape and mix the contents by moving the strip holder in a circular motion on the benchtop for 60 seconds. Be careful not to spill the contents.
8. Incubate 30 minutes at room temperature.
9. Remove the covering and decant the contents of the wells into a sink. Wash the strips **three times** using the **diluted washing buffer solution**. Please use at least a volume of 250 µL of washing buffer for each well in each washing step. Remaining buffer in the wells should be removed by patting the inverted plate dry on a stack of paper towels.
10. Add 100 µL of **substrate (color) solution** to the wells. Cover the wells with parafilm or tape and mix the contents by moving the strip holder in a circular motion on the benchtop for 60 seconds. Be careful not to spill the contents. Incubate the strips for 20 minutes at room temperature. Protect the strips from direct sunlight.
11. Add 50 µL of **stop solution** to the wells in the same sequence as for the substrate solution.
12. Read the absorbance at 450 nm using a microplate ELISA photometer within 15 minutes after the addition of stopping solution.

G. Evaluation

The evaluation of the ELISA can be performed by comparing the raw absorbance value for each control and sample, averaged between the two (or three) replicates, to the following known values. For both dilutions, raw absorbance (averaged between replicates) values of the kit positive control should range from 0.888 to 1.332; values for the kit negative control should range from 0.108 to 0.162. Absorbance values outside this range may indicate an invalid test. Uninoculated media, incubated without dilution alongside samples per USDA guidelines, should return raw absorbance values (averaged between replicates) ranging from 0.147 to 0.220. Values outside this range may indicate contamination.

Screening dilution: for samples incubated using the 1:4 dilution, absorbance values higher than 0.205 should be considered positive and submitted to further testing. Sensitivity using this method is as low as 30 pg/mL (approximately n cells per 25 g beef).

Outbreak Related dilution: for samples incubated using the 1:10 dilution, absorbance values higher than 0.215 should be considered positive and submitted to further testing. Sensitivity using this method is as low as 30 pg/mL (approximately n cells per 10 g beef).



Safety Data Sheet

Section 1: Product and Company Identification

1.1 Product Identifiers:

Product Name: Shiga Toxin 1 (STX 1), Shiga Toxin 2 (STX 2) ELISA Plate Kits

Product Code: 542000, 542010

1.2 Identified Use: Determination of Shiga Toxin 1 and 2 in samples. **Restrictions on Use:** For research use only.

1.3 Company: Abraxis, Inc., 124 Railroad Drive, Warminster, PA 18974 USA, info@abraxiskits.com +1(215) 357-3911, FAX +1(215) 357-5232

1.4 Emergency Telephone Number: +1(215) 357-3911

Section 2: Hazard(s) Identification

2.1 Classification of the mixture: Not a hazardous mixture.

2.2 GHS Label elements, including precautionary statements: Not applicable.

2.3 Hazards not otherwise classified (HNOC) or not covered by GHS: None known.

2.4 Unknown acute toxicity: None known.

Section 3: Composition / Information on Ingredients

3.2 Mixtures: *Contains no hazardous ingredients at levels requiring disclosure by the OSHA Hazard Communication Standard (29 CFR 1910.1200), however it contains minor amounts of materials considered hazardous. We recommend handling all substances with caution.*

Section 4: First Aid Measures

4.1 Description of first aid measures: Consult a physician. Show this safety data sheet to the doctor in attendance. Move out of dangerous area.

If inhaled: If breathed in, move person into fresh air. If not breathing, give artificial respiration. Consult a physician.

In case of skin contact: Wash off with soap and plenty of water. Consult a physician.

In case of eye contact: Rinse thoroughly with plenty of water for at least 15 minutes and consult a physician.

If swallowed: Never give anything by mouth to an unconscious person. Rinse mouth with water. Consult a physician.

4.2 Most important symptoms and effects, both acute and delayed: No data available

4.3 Indication of any immediate medical attention and special treatment needed: No data available. Treat symptomatically.

Section 5: Fire-fighting Measures

5.1 Suitable extinguishing media: Use an extinguishing agent suitable for the surrounding fire.

5.2 Special hazards arising from the substance or mixture: None known

5.3 Advice for firefighters: Wear self-contained breathing apparatus for fire-fighting if necessary.

5.4 Further information: No data available

Section 6: Accidental Release Measures

6.1 Personal precautions, protective equipment and emergency procedures: Use personal protective equipment (see section 8). Avoid dust formation. Avoid breathing vapors, mist, dust, or gas. Ensure adequate ventilation. Evacuate personnel to safe areas.

6.2 Environmental precautions: Prevent further leakage or spillage if safe to do so. Do not let product enter drains. Discharge into the environment must be avoided.

6.3 Methods and materials for containment and cleaning up: Solids (if applicable): Pick up and arrange disposal without creating dust. Sweep up and shovel. Liquids (if applicable): Absorb with liquid-binding material (sand, diatomite, acid binders, universal binders, sawdust). Keep in suitable, closed containers for disposal.

6.4 Reference to other sections: For information on safe handling see section 7.

For information on personal protection see section 8.

For information on disposal see section 13.

Section 7: Handling and Storage

7.1 Precautions for safe handling: See section 2. Avoid inhalation of vapors and contact with skin and eyes. Wear appropriate personal protective equipment. Do not eat, drink, or smoke in work area.

7.2 Precautions for safe storage: Keep container(s) tightly closed in a dry, well-ventilated place. Protect from physical damage. See label or product insert for appropriate storage temperature and additional specific information.

7.3 Specific end use(s): No data available

Section 8: Exposure Controls / Personal Protection

8.1 Control parameters: Not applicable.

8.2 Exposure controls:

Appropriate engineering controls: Provide adequate ventilation. Handle in accordance with good industrial hygiene and safety practice. Wash hands before breaks and at the end of workday. Keep away from food and beverages.

Personal protective equipment: The usual precautionary measures, including eye/face/skin protection, should be taken when handling any chemical. Avoid contact with eyes, skin, and clothing.

Eye protection: As with handling of any chemical, wear approved safety goggles.

Skin protection: Handle with gloves. No specific information regarding glove material or thickness is available, but gloves must be impermeable and resistant to the substance being handled. Gloves must be inspected prior to use. Use proper glove removal technique (without touching glove's outer surface) to avoid skin contact with this product. Dispose of contaminated gloves after use in accordance with applicable laws and good laboratory practices. Wash and dry hands.

Respiratory protection: As with any chemical, where excessive vapor, mist, or dust may result, use a chemical fume hood or approved respiratory protection equipment.

Body protection: Lightweight, protective clothing.

Section 9: Physical and Chemical Properties

9.1 Information on basic physical and chemical properties of the mixture

Appearance: Multiple

Odor: Characteristic

Odor Threshold: No data available

pH: Multiple

Melting point/freezing point: No data available

Initial boiling point and boiling range: No data available

Flash point: Not applicable

Evaporation rate: No data available

Flammability (solid, gas): No data available

Upper/lower flammability or explosive limits: No data available

Vapor pressure: No data available

Vapor density: No data available

Relative density: No data available

Water solubility: Various

Partition coefficient: n-octanol/water: No data available

Auto-ignition temperature: Not applicable

Decomposition temperature: No data available

Viscosity: No data available

Explosive properties: No data available

Oxidizing properties: No data available

9.2 Other information: No data available

Section 10: Stability and Reactivity

10.1 Reactivity: No data available

10.2 Chemical stability: Stable under recommended storage conditions.

10.3 Possibility of hazardous reactions: No data available

10.4 Conditions to avoid: No data available

10.5 Incompatible materials: No data available

10.6 Hazardous decomposition products: No data available. In the event of fire: see section 5.

Section 11: Toxicological Information

11.1 Information on toxicological effects

Acute toxicity: Not available. To the best of our knowledge, the chemical, physical, and toxicological properties of this product have not been thoroughly investigated.

Inhalation: No data available **Ingestion:** No data available

Skin contact: Irritant to skin and mucous membranes.

Eye contact: May cause eye irritation in susceptible persons.

Respiratory or skin sensitization: No data available

Aspiration hazard: No data available

Mutagenicity: No data available

Carcinogenicity

IARC: No component of this product present at levels greater than or equal to 0.1% is identified as probable, possible or confirmed human carcinogen by IARC.

ACGIH: No component of this product present at levels greater than or equal to 0.1% is identified as a carcinogen or potential carcinogen by ACGIH.

NTP: No component of this product present at levels greater than or equal to 0.1% is identified as a known or anticipated carcinogen by NTP.

OSHA: No component of this product present at levels greater than or equal to 0.1% is identified as a carcinogen or potential carcinogen by OSHA.

Teratogenicity: No data available

Reproductive/fertility toxicity: No data available

Specific target organ toxicity, single exposure: No data available

Specific target organ toxicity, repeated exposure: No data available

Section 12: Ecological Information

12.1 Toxicity: No data available

12.2 Persistence and degradability: No data available

12.3 Bioaccumulative potential: No data available

12.4 Mobility in soil: No data available

12.5 Results of PBT and vPvB assessment: No data available

12.6 Other adverse effects: An environmental hazard cannot be excluded in the event of unprofessional handling or disposal.

Section 13: Disposal Considerations

13.1 Waste treatment methods

Product: All waste must be handled and disposed according to local, state, and federal regulations. Avoid disposing large volumes in sewer.

Contaminated packaging: All waste must be handled and disposed according to local, state, and federal regulations.

Refer to sections 7 and 8 for safe handling guidance.

Section 14: Transport Information

UN Number: Not regulated

UN Proper shipping name: Not classified as dangerous in the meaning of transport regulations.

Transport hazard class(es): No data available

Packing group: No data available

Environmental hazard: No data available

Bulk transport: No data available

Special considerations: No data available

Section 15: Regulatory Information

To the best of our knowledge, this product contains no substances which, at their given concentrations, are considered hazardous by other regulatory agencies. Refer to section 3.

Section 16: Other information

This information is based on our present knowledge. While Abraxis, Inc. believes that the data contained herein are factual and the opinions expressed represent a best effort to present accurate information, the data are not to be taken as a warranty or representation for which Abraxis, Inc. assumes legal responsibility. The information shall not be taken as being all-inclusive and is to be used only as a guide. The data are offered solely for the user's consideration, investigation, and verification. These suggestions should not be confused with either state, municipal, or insurance requirements, or with national safety codes and constitute no warranty. Any use of these data and information must be determined by the user to be in accordance with applicable federal, state, and local regulations.

All materials and mixtures may present unknown hazards and should be used with caution. Since Abraxis, Inc. cannot control the methods, volumes, or conditions of use of this product, Abraxis, Inc. shall not be held liable for any damages or losses resulting from the handling or from contact with the product as described herein. An individual technically qualified to handle potentially hazardous chemicals must supervise the use of this material. This product is sold for research use only. It is not for any human or animal therapeutic or clinical diagnostic use.

Date this SDS was prepared: 5/24/2016

Version: 2

Changes from previous version: Abraxis, LLC changed to Abraxis, Inc.